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The TS-100C Pro is a programmable laboratory thermo-shaker that combines precise heating and cooling, intensive mixing and interchangeable thermoblocks in one compact instrument. Designed for repeatable sample-preparation workflows, it allows users to create and save multi-step programs directly on the touchscreen. This reduces repeated manual parameter changes and the need to transfer samples between separate heating, cooling and mixing instruments.

KEY BENEFITS

- **Heating, cooling and mixing** in one instrument for efficient temperature-controlled sample preparation.
- **Mixing speeds up to 2500 rpm** support intensive mixing, bead resuspension, extraction and lysis workflows.
- **Automatic thermoblock recognition** identifies the installed block and limits the maximum mixing speed accordingly.
- **Quick thermoblock exchange** without external tools or separate cable connections reduces setup time between applications.
- **Advanced on-screen programming** with up to 20 saved programs and up to 20 steps per program, including holds, repetitions and parameter transitions.
- **Clear touchscreen interface** displays set and actual temperature, speed, time and program status at a glance.
- **Quick Mix function** starts a short mixing cycle using the selected parameters without creating a complete program.
- **Optional heated lid** helps minimise condensation during temperature-controlled workflows. The instrument recognises the lid automatically, while lid heating can be switched on or off as required.
- **Five configurable user profiles and screen lock** help protect established methods and prevent unintended changes in shared laboratories.
- **Integrated system and maintenance information** provides access to the serial number, firmware version, operating hours, calibration status, service reminders and diagnostics. The operating manual can be accessed using an on-screen QR code.

APPLICATIONS

The TS-100C Pro is suitable for **DNA and RNA extraction, NGS library preparation, magnetic-bead cleanup and resuspension, cell lysis, protein extraction, enzymatic reactions, incubation, hybridisation and analytical sample preparation.**

A range of interchangeable thermoblocks allows the instrument to be configured for different compatible tubes, strips and plate formats.

*Maximum mixing speed depends on the installed thermoblock.

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