

FTA-1 & FTA-2i

Aspirators with trap flask



Contents

1.	About this edition of user instructions.....	3
2.	Safety precautions	4
3.	General information.....	5
4.	Getting started.....	6
5.	Operation	9
6.	Specifications	10
7.	Care and maintenance	12
8.	Warranty.....	14
9.	EU Declaration of conformity.....	15

1. About this edition of user instructions

The current edition of the user manual applies to the following models and versions of aspirators with trap flasks:

- **FTA-1** version V.4AW
- **FTA-2i** version V.1A01

2. Safety precautions



Caution!

Make sure you have fully read and understood the present Manual before using the equipment. Please pay special attention to sections marked by this symbol.

GENERAL SAFETY

- The protection provided can be ineffective if the operation of the appliance does not comply with the manufacturer's requirements.
- Save the unit from shocks and falling.
- Store and transport the unit at ambient temperatures between -20°C and +60°C and maximum relative humidity of 80%.
- After transportation or storage and before connecting it to the electric circuit, keep the unit under room temperature for 2-3 hrs.
- Use only original parts and accessories, provided by manufacturer for this product.
- Before using any cleaning or decontamination methods except those recommended by the manufacturer, check with the manufacturer that the proposed method will not damage the equipment.
- Do not make modifications in design of the unit.

ELECTRICAL SAFETY

- Connect only to the mains with voltage corresponding to that on the serial number label.
- Use only the external power supply provided with this product.
- Ensure that the power plug is easily accessible during use.
- Disconnect the unit from the mains before moving.
- If liquid penetrates into the unit, disconnect it from the mains and have it checked by a repair and maintenance technician.
- Do not operate the unit in premises where condensation can form. Operating conditions of the unit are defined in the **Specifications** section.

DURING OPERATION

- Do not operate the unit in environments with aggressive or explosive chemical mixtures. Please contact manufacturer for possible operation of the unit in specific atmospheres.
- Do not operate the unit if it is faulty or has been installed incorrectly.
- Do not use outside laboratory rooms.
- Do not use the unit for aggressive organic compounds, chlorinated solvents and strong conjugated acids.

BIOLOGICAL SAFETY

- It is the user's responsibility to carry out appropriate decontamination if hazardous material is spilt on or penetrates into the equipment.
- Empty the trapping flask when the liquid level reaches maximum. Take necessary precautions utilizing waste liquid in accordance with general laboratory standards.

3. General information

FTA-1 & FTA-2i aspirators with trap flask have been designed for liquid removal from micro-test tubes (e.g. for supernatant removal during DNA/RNA extraction), ELISA plates (e.g. performing several repeated well washing cycles), cell culture flasks (cell media exchange and washing during cell growth) and other vessels during various laboratory techniques.

Model **FTA-1** has a 1 L glass trap flask and is designed to be a low-cost solution. Aspiration is possible with a single standard 200 µL tip and the 8-tipped **MA-8** adapter.

Model **FTA-2i** comes with a 2 L polypropylene trap flask and a liquid overflow sensor that stops collected material from damaging the filter. Aspiration is possible with single 200 and 1000 µL tip using the **MA-U** adapter or with the handheld vacuum controller **HAS-1** with different attachments such as 8-channel adapter with stainless steel tips, 8-channel adapter for 200 µl tips, single channel adapter for 200 µl tips and others.

Microbiologic suction filter eliminates risks of contamination caused by bacteria, viruses, and infected particles from samples. Such contamination can be distributed by suction pump or central vacuum units. Microbiological suction filter is hydrophobic and has very high contamination removal bacterial efficiency – up to 99.99% of particles bigger than 0.027 micrometres (e.g. Hepatitis A, B and C, HIV, etc.).

In order to prolong performance and efficiency of device it is not allowed to use aggressive organic compounds (e.g. phenol) and strong conjugated acids (e.g. guanidine) for aspiration. Please, contact manufacturer in case of any questions regarding operation of device.

4. Getting started

- 4.1. **Unpacking.** Remove packing materials carefully and retain them for future shipment or storage of the unit. Examine the unit carefully for any damage incurred during transit. The warranty does not cover in-transit damage. Warranty covers only the units transported in the original package.
- 4.2. **Complete set.** Package contents:
- 4.2.1. **FTA-1:**
- FTA-1, aspirator with trap flask 1 pce.
 - Hydrophobic filter 1 pce.
 - Aspiration tips, 200 µL 50 pcs.
 - Holder tubes (1.5 ml) 10 pcs.
 - Spare hydrophobic filter 1 pce.
 - External power supply 1 pce.
 - Operating manual, declaration of conformity 1 copy
 - **MA-8**, 8-channel aspiration tip with holderon request
- 4.2.2. **FTA-2i:**
- FTA-2i, aspirator with flask trap 1 pce.
 - **MA-U**, universal adapter 1 pce.
 - Hydrophobic filter 1 pce.
 - Aspiration tips, 200 µL 50 pcs.
 - Aspiration tips, 1000 µL 10 pcs.
 - Holder tubes (2 ml) 5 pcs.
 - Spare hydrophobic filter 1 pce.
 - External power supply 1 pce.
 - Operating manual, declaration of conformity 1 copy
 - **MA-8**, 8-channel aspiration tip with holderon request
 - **HAS-1**, hand operator set with SH-6 standon request

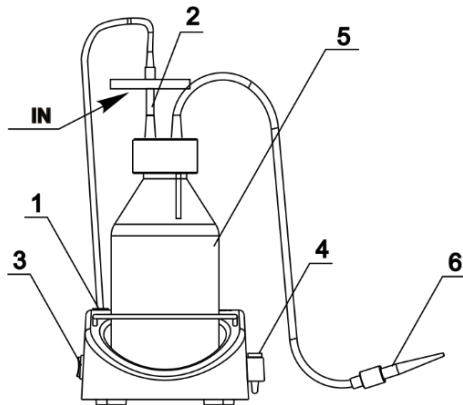


Figure 1. Assembled FTA-1

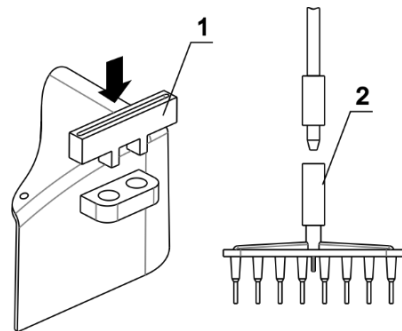


Figure 2. Installation of MA-8 aspiration tip

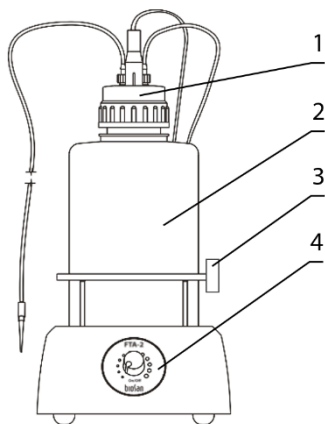


Figure 3. FTA-2i front view

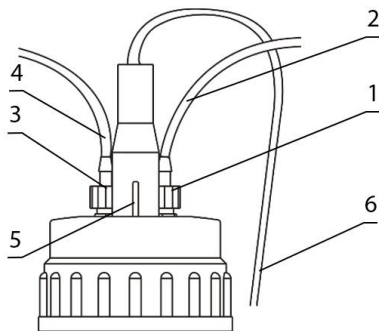


Figure 4. FTA-2i lid with sensor and tubing



Figure 5. MA-U, universal adapter

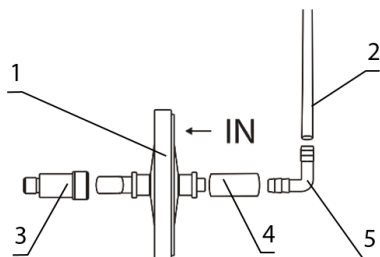


Figure 6. FTA-2i filter setup

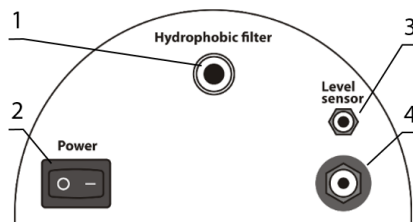


Figure 7. FTA-2i rear panel

4.3. FTA-1 setup.

- Place the unit on horizontal even working surface.
- Install the trapping flask (fig. 1/5) on the unit.
- Assemble the lid, aspiration tubing and filter set (fig. 1/2). Attach aspiration tip (fig. 1/6) to the tubing.



Caution! Position the filter with the **IN** marking pointing towards the unit (fig. 1/2).

- Screw on the lid with the filter set and aspiration tubing to trap flask. Ensure that the filter set is tightly inserted.
- Connect the filter tubing to the pump (fig. 1/1) located at the top of the base behind the trap flask, by turning the fitting clockwise.
- Place the microtubes for aspiration tip into the holder (fig. 1/4).
- Connect the external power supply unit into the socket at the rear side of the unit and position the unit for an easy access to the external power supply and the power switch.

- 4.3.1. **Optional MA-8 aspiration tip installation.** Remove microtubes and aspiration tip from the holder (fig. 1/4). Place MA-8 storing holder (fig. 2/1) into the holder. Remove aspiration tip (fig. 1/6) from the aspiration tubing and connect the **MA-8** aspiration tip (fig. 2/2). Store the aspiration tip in the new holder.

4.4. FTA-2i setup.

- Place the unit on horizontal even working surface.
- Install the trapping flask (fig. 3/2) on the unit.
- Screw on the lid with the sensor needles and valves for tubing (fig. 3/1) to trapping flask.
- Assemble the filter set (figure 6) and connect it to the port on the rear panel of the unit (fig. 7/1).



Caution! Position the filter with the **IN** marking pointing away from the unit (fig. 6/1).

- Attach the fitting (fig. 4/1) of the tubing to the filter (fig. 4/2) to the valve with **white** ring on the lid.
- Ensure that the filter set is tightly fixed in the slot (fig. 7/1).
- Assemble the **MA-U** aspiration tip set (figure 5).
- Attach the fitting (fig. 4/3) of the tubing for liquids (fig. 4/4 and 5/3) to the valve with the **blue** ring on the lid.
- Connect the sensor cable (fig. 4/5) to the lid of the unit and to the port on the rear panel of the unit (fig. 5/3).
- Place the microtubes for aspiration tip into the holder (fig. 3/3).
- Connect the external power supply unit into the socket at the rear side of the unit (fig. 7/4) and position the unit for an easy access to the external power supply and the power switch.

4.4.1. **Optional MA-8 aspiration tip installation.** Remove microtubes and aspiration tip from the holder (fig. 3/3). Remove the aspiration tip (fig. 5/1) from the **MA-U** adapter and connect the **MA-8** aspiration tip (fig. 8/1).

4.4.2. **Optional HAS-1 hand operator installation.** Unscrew the **MA-U** adapter from the aspiration tube (fig. 5/3) and connect the **HAS-1** hand operator (fig. 8/6) by screwing it to the fitting.

5. Operation

- 5.1. Connect the external power supply to the mains.
- 5.2. Turn the **Power** switch (fig. 1/3 or 7/2) to the position **I** (on). The pump starts working.
- 5.3. (For model **FTA-2i**) Regulate the aspiration intensity with the regulator knob on the front panel of the unit (fig. 3/4). Clockwise turning direction increases power, counter clockwise – decreases. Current intensity is indicated by a light ring around the regulator knob.
- 5.4. Touch the surface of the liquid with the aspiration tip (fig. 1/6, 5/1 or **MA-8**) to remove the liquid.



Caution!

During operation, do not allow the collected liquid to exceed the maximum level. Empty the trapping flask when the liquid level reaches maximum allowed level, according to **5.7**.



Note.

In model **FTA-2i**, when collected liquid exceeds the maximum level, the sensor turns off the pump and produces sound and light warning signals.

- 5.5. (For model **FTA-2i**) When using the optional **HAS-1** hand regulator, touch the liquid with the aspiration tip, then press and hold the blue button on the regulator to collect.
- 5.6. After finishing the operation, switch off the **Power** switch and disconnect the external power supply from the mains.
- 5.7. **Emptying the trapping flask.** Turn off the unit. Remove the tubings and cables from the lid. Remove the closed flask from its place. Unscrew the lid and empty the trapping flask. Screw the lid back on and return the trapping flask on the unit. Replace the tubings and the cables.



Caution!

Take necessary precautions in utilizing waste liquid in accordance with general laboratory standards. For information regarding autoclaving procedures, see **7.4**.

6. Specifications

The unit is designed for operation in cold rooms, incubators (except CO₂ incubators) and closed laboratory rooms at ambient temperature from +4°C to +40°C in a non-condensing atmosphere and maximum relative humidity 80% for temperatures up to 31°C decreasing linearly to 50% relative humidity at 40°C.

Biosan is committed to a continuous programme of improvement and reserves the right to alter design and specifications of the equipment without additional notice.

	FTA-1	FTA-2i
Aspiration speed	up to 1.7 mL/s	up to 27.8 mL/s
Vacuum regulation	None	Yes
Working vacuum pressure	-500 mbar	-200 ... -800 mbar
Overflow protection sensor	None	Yes, invasive
Overflow protection actions	None	Pump stop, sound and light signals
Trapping flask	1 L, glass	2 L, polypropylene (PP)
Air filtration	Hydrophobic microbiologic air filter 2200/02 ¹	
Filter pore diameter	0.027 µm	
Input current	DC 12 V, 300 mA	DC 12 V, 1 A
Power consumption	3.6 W	10.8 W
External power supply	input AC 100-240 V, 50/60 Hz; output DC 12 V	
Dimensions	160x210x340 mm	185x290x390 mm
Weight ²	1.7 kg	1.85 kg

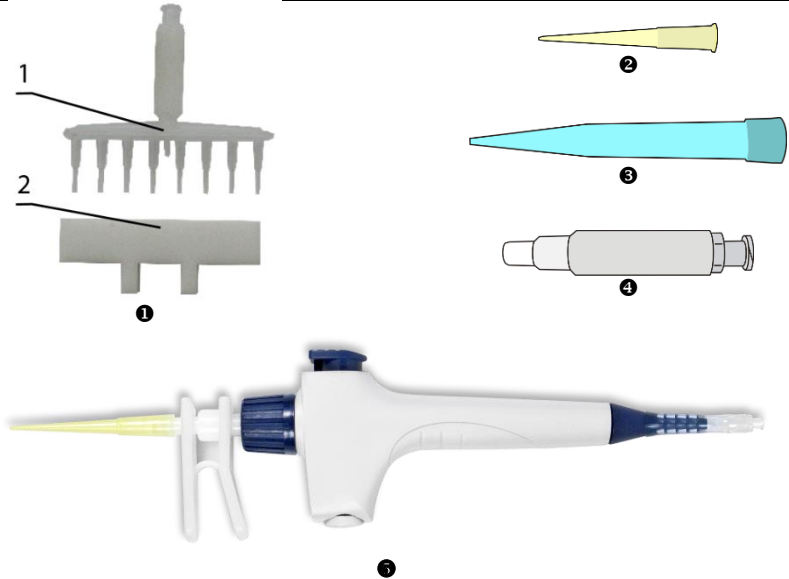


Figure 8. Optional accessories and replacement parts

¹ Eliminates risk of contamination from the trapping flask with bacteria, viruses and infected particles
² Accurate within ±10 %

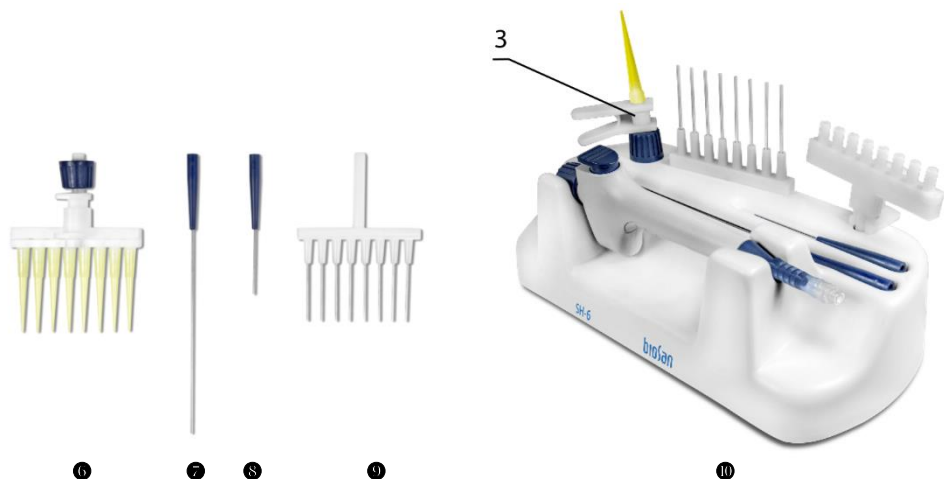


Figure 8. Optional accessories and replacement parts (continued)

Table 1. Optional accessories

Description	Catalogue number
MA-8 , 8-channel aspiration tip with holder (fig. 8/1 1)	BS-040108-BK
HAS-1 , hand operator set (only for FTA-2i). Set includes: Handheld vacuum controller with adapter (fig. 6/5), 8-channel adapter (with ejector) for 200 μ L tips (fig. 6/6), 1-channel adapter with 125 mm stainless steel pin (fig. 6/7), 1-channel adapter with 40 mm stainless steel pin (fig. 6/8), 8-channel adapter with 35 mm stainless steel pin (fig. 6/9), 1-channel adapter (with ejector) for 200 μ L tips (fig. 6/3), Stand SH-6 for hand operator parts (fig. 6/10)	BS-040118-PK

Table 2. Replacement parts for FTA-1

Description	Catalogue number
Hydrophobic filter	BS-040108-S25
Aspiration tip, 200 μ L, 1000 pcs. (fig. 8/2)	BS-000001-S04
Aspiration tip, 8 channels, for MA-8 (fig. 8/1)	BS-040108-S24
Trapping flask, 1 L, glass, with lid, inserts and sealing ring	BS-040108-S11
Tubing and fitting set	BS-040108-S22

Table 3. Replacement parts for FTA-2i

Description	Catalogue number
MA-U , universal adapter for 200/1000 μ L single use tips (fig. 8/4)	BS-040118-AK
Hydrophobic filter	BS-040108-S25
Aspiration tip, 200 μ L, 1000 pcs. (fig. 8/2)	BS-000001-S04
Aspiration tip, 1000 μ L, 1000 pcs. (fig. 8/3)	BS-000001-S05
Aspiration tip, 8 channels, for MA-8 (fig. 8/1)	BS-040108-S24
Trapping flask, 2 L, polypropylene, with lid, inserts and sealing ring	BS-040120-S11
Tubing and fitting set	BS-040120-S14
Sensor cable	BS-040120-S12

7. Care and maintenance

7.1. If the unit requires maintenance, disconnect the unit from the mains and contact Biosan or your local Biosan representative.

7.2. All maintenance and repair operations must be performed only by qualified and specially trained personnel.

7.3. Cleaning and disinfection.

7.3.1. Glass and plastic outer parts. Standard ethanol (75%) or other cleaning agents recommended for cleaning of laboratory equipment can be used for cleaning and decontamination.

7.3.2. Silicone and internal parts. Use only distilled water, no less than 0.5 L.

7.4. Autoclaving procedures.

	Lid, fittings, tubing	Flask	Hydrophobic filter, cables, base pump
FTA-1	121°C and 1 atm., 15 min.	121°C and 1 atm., 55 min.	Do not autoclave
FTA-2i		121°C and 1 atm., 15 min.	



Caution!

Do not put a closed flask in the autoclave, as it might irreversibly deform. Completely disengage the thread on the cap.



Note.

Hydrophobic filter and sensor cable are sensitive to moisture and are not autoclavable.

Tubing is made of silicone and deteriorate after frequent autoclaving and should be replaced.

7.5. **Tubing replacement.** Change the silicone tubing every 6 months.

7.6. **Filter replacement.** Replace the filter as soon as it accumulates visible traces of dirt, or when the liquid is sucked into it. Maximum working time for the filter is 30 days. When working with infected materials, replace the filter immediately after finishing the procedure to prevent the spread of infection.

7.6.1. Changing procedure. Remove the filter (fig. 1/2 or 6/1) by detaching it from adapters. Insert a new filter and reconnect the adapters.



Caution!

Observe the orientation of the filter:

- For model **FTA-1**, the **IN** marking should be facing towards the unit (figure 1).
- For model **FTA-2i**, the **IN** marking should be facing away from the unit (figure 6).

7.7. Troubleshooting.

Problem	Possible reason	Troubleshooting
Unit not working	No power	Check external power supply and connector
	(FTA-2i) Trapping flask full	Sensor shut down the motor. Empty the trapping flask according to 5.7.
	(FTA-2i) Sensor not connected properly	Check sensor cable and attachments
Aspiration slow or absent	(FTA-2i) Aspiration intensity set to small	Turn the regulator knob clockwise to increase aspiration.
	Breach in the system	Check all the tubing and fittings. Tightly close the lid of the unit
	Dirt in the tubing	Clean the tubing, see 7.3.
	Liquid in the filter	Replace the filter
(FTA-2i) Sensor activated, but the flask is not full	Foam in the trapping flask activated the sensor	Do not use the unit with a foaming liquid or use antifoaming agent.

8. Warranty

- 8.1. The Manufacturer guarantees the compliance of the unit with the requirements of Specifications, provided the Customer follows the operation, storage and transportation instructions.
- 8.2. The warranted service life of the unit from the date of its delivery to the Customer is 24 months. For extended warranty, see **8.5**.
- 8.3. Warranty covers only the units transported in the original package.
- 8.4. If any manufacturing defects are discovered by the Customer, an unsatisfactory equipment report shall be compiled, certified and sent to the local distributor address. To obtain the claim form, visit section **Technical support** on our website at link below.
- 8.5. Extended warranty.
- For **FTA-2i**, the *Premium* class model, one year of extended warranty is available free of charge after registration, during 6 months from the date of sale. Online registration form can be found in section **Warranty registration** on our website at the link below.
 - For **FTA-1**, the *Basic Plus* class model, extended warranty is a paid service. Contact your local Biosan representative or our service department through the **Technical support** section on our website at the link below.
- 8.6. Description of the classes of our products is available in the **Product class description** section on our website at the link below.

Technical support



biosan.lv/en/support

Registration



biosan.lv/register-en

Product class description



biosan.lv/classes-en

- 8.7. The following information will be required in the event that warranty or post-warranty service comes necessary. Complete the table below and retain for your records.

Model	FTA-1 & FTA-2i , aspirators with trap flask
Serial number	
Date of sale	

9. EU Declaration of conformity

EU Declaration of Conformity

Unit type Aspirators with trap flask

Models **FTA-1, FTA-2i**

Serial number 14 digits styled XXXXXYYMMZZZZ, where XXXXXX is model code, YY and MM – year and month of production, ZZZZ – unit number.

Manufacturer SIA BIOSAN
Latvia, LV-1067, Riga, Ratsupites str. 7/2

Applicable Directives EMC Directive 2014/30/EU
LVD Directive 2014/35/EU
RoHS2 2011/65/EU
WEEE 2012/19/EU

Applicable Standards LVS EN 61326-1: 2013
Electrical equipment for measurement, control and laboratory use. EMC requirements. General requirements.
LVS EN 61010-1: 2011
Safety requirements for electrical equipment for measurement, control, and laboratory use. General requirements.

We declare that this product conforms to the requirements of the above Directives




Signature

Svetlana Bankovska
Managing director

19.07.2016.

Date

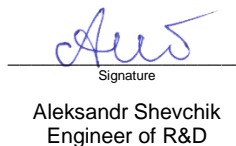

Signature

Aleksandr Shevchik
Engineer of R&D

19.07.2016

Date

Biosan SIA

Ratsupites 7, build.2, Riga, LV-1067, Latvia
Phone: +371 67426137 Fax: +371 67428101
<http://www.biosan.lv>

Edition 1.-4.01 – March 2018